### **EQUIPMENT**

1.	TEST PLATES: smooth polished surface stainless steel or aluminum test plates (records)
	a) plate diameter 8.00 ± 0.02 mm or 25.00 ± 0.05 mm dia
	i) with raised portion of minimum 1.50 mm
	b) top and bottom plates have same diameter with minimum thickness ≥ 1.50 mm
2.	ENVIRONMENTAL CHAMBER:
	a) controlled temperature test chamber (by liquid or gas to 0.1°C)
	b) if cooled by air of gas a suitable drier is used to prevent condensation or ice formation
	c) if cooled by liquid circulating bath used
	d) chamber completely encloses the top and bottom plates
3.	TEMPERATURE CONTROLLER: maintains specimen temperature from 3 to 88°C
	within 0.1°C
4	INTERNAL TEMPERATURE DETECTOR: platinum resistance thermometer (PRT) in
	intimate contact with fixed plate, controls temperature in range of 3 to 88°C to 0.1°C and
_	provides continuous readout <i>(records)</i>
5	a) reference thermometer NIST traceable liquid in glass or NIST traceable electronic
	thermometric device and maintained as the laboratory standard — shall be used to
	standardize portable thermometer (records)
	i) liquid-in-glass within an ice point and standardized according to ASTM E563
	ii) portable thermometer having a with accuracy of ± 0.05°C and resolution of 0.01°C,
	calibrated at least yearly using NIST-traceable reference standard in accordance with
	ASTM E77
	b) electronic thermometer with resistive detector and digital readout (may be the reference
	thermometer if it fits within dummy specimen) calibrated at lest yearly (records)
6.	LOADING DEVICE: applies sinusoidal oscillatory load at frequency of 10.0 ± 0.1 rad/sec
	a) if strain controlled, shall apply a cyclic torque sufficient to cause an angular rotation
	strain to within 100 µ rad of the strain specified
	b) if stress controlled, shall apply a cyclic torque accurate to within 10 mN.m of the torque
	specified
6	TEMPERATURE CORRECTION: must be traceable to NIST standards (records)
0.	a) silicone wafer of thickness or dia. < 2.0 mm used for dummy specimen
	b) dummy test specimen used
7.	DATA ACQUISITION SYSTEM: (demo)
	a) provides a record of
	i) temperature to 0.1°C
	ii) frequency to 1 %
	iii) torque to 10 mN.m
	iv)deflection angle to 100 µrad
	b) measures and records
	i) complex shear modulus (G*) in range of 100Pa to 10MPa with accuracy of 1.0%
	or less

	ii) phase angle (D) in range of 0 to 90° (with accuracy of 0.1°)
	iii) calculates shear stress, shear strain, complex shear modulus and phase angle
9.	SPECIMEN MOLD: silicone mold of defined dimensions (optional)
	.CLEANING SOLVENTS: mineral oil, toluene, citrus-based, mineral spirits; acetone to
	remove solvent residue
	.WIPING MATERIAL: clean cloth, paper towel, cotton swabs
12	. 3F LOIMEN TRIMINER. Metal spatula of similar with straight edge 4 mm wide
	VERIFICATION AND CALIBRATION
	TEMPERATURE: dummy asphalt binder prepared or silicone wafer used
2.	TEMPERATURE CORRECTION: (demo – records)
	a) temperature measured using dummy sample or silicone wafer with the reference thermal detector and compared to DSR RTD temperature
	b) temperature at equilibrium (difference between reference detector and DSR RDT
	over a five min. period)
	c) reference detector temperature vs DSR temperature plotted at 6 °C intervals over
	range of test temperatures
	<ul> <li>d) appropriate correction applied to DSR RDT if not in agreement within ± 0.1°C</li> <li>i) does softwear incorporate temperature corrections</li> </ul>
3	DSR TORQUE TRANSDUCER:using
٥.	a) standard viscosity with DSR complex viscosity (G*/(rad/s) within 3.0 % of capillary
	viscosity of the liquid standard
	b) using supplied fixtures if available
4.	ANGULAR DISPLACEMENT TRANSDUCER: calibrated as per directions and fixtures
	supplied with apparatus (if available – otherwise not required)
5.	TEST PLATE DIAMETERS:
	a) 8.00 ± 0.02 mm and 25.0 ± 0.05
	b) if in range of 7.9 to 8.1 mm and 24.5 to 25.5 mm diameters measured and maintained in log
6	OVERALL CALIBRATION: verified using reference viscosity fluids (demo – records)
	VERIFICATION: is carried out every 6 months or when DSR is
•	a) newly installed
	b) relocated
	c) accuracy suspected
	DDEDADATION OF ADDADATHO (Acces)
4	PREPARATION OF APPARATUS (demo)
	APPARATUS: set up as per manufacturer recommendation  TEST PLATES: checked for damage, cleaned and mounted on test fixtures and
۷.	tightened firmly
3.	TEST TEMPERATURE: selected according to grade of asphalt binder(M320) and
	equipment temperature stabilized to within 0.1 °C of target temperature
	Page 2 of 5

4.	GAP SETTINGS:
	<ul><li>a) zero gap set (at temperature) by:</li><li>i) manually by spinning moveable plate while closing the gap until the moveable plate touches the fixed plate</li></ul>
	ii) automatically closing the gap while watching the normal force, after contact set zero gap approx. to zero force
	b) plates moved apart to approximate test gap settings
	c) plates preheated to:
	i) test temperature for 25 mm plates (or lowest temperature if testing at more than one temperature)
	ii) to 34 to 46°C range for 8 mm plates
	II) to 54 to 40 o range for 6 min plates
	PREPARATION OF SAMPLES AND TEST SPECIMEN (demo)
1.	TEST SPECIMENS:
	a) binder heated until just fluid enough to be transferred (> 135°C to be avoided
	if possible)
	i) to removable plate removed, sample poured continuously approximately 15 mm
	above plate until plate covered to within approximately 2 mm of the perimeter, allowed to stiffen and plate placed in the DSR
	ii) to one of the plates using a rod or spatula
	iii) by pouring the binder into a silicone rubber mold, cooled, demolded and centered
	on plate
	b) immediately after samples placed on the plates, the plates are moved together to:
	i) 1 mm plus 0.05 mm for 25 mm plate
	ii) 2 mm plus 0.10 mm for 8 mm plate
	c) gap decreased to produce bulge, typically: i) 0,05 mm for 25 mm plate
	ii) 0.10 mm for 8 mm plate
	d) excess binder trimmed flush to plates
	,
	<u>PROCEDURE</u>
1.	TEMPERATURE:
	a) test temperature set to ± 0.1°C, including any offset
	b) desired temperature (including any offset) is maintained at the test temperature ± 0.1°C for at least 10 min before start
2	STRAIN CONTROL MODE: controlled to within 20 % of target value as per table below
	STRESS CONTROL MODE: controlled to within 20 % of target value as per table below

Page 3 of 5

	WHEN TESTING FOR COMPLIANCE WITH M320G						
Ta	arget strain Value	es	Target Stress Values				
Material	Target, %	Range, %	Material	Target, %	Range, %		
Original	12	9 to 15	Original	0.12	0.09 to 0.15		
RTFO	10	8 to 12	RTFO	0.22	0.18 to 0.26		
Residue			Residue				
PAV Residue	1	0.8 to 1.2	PAV Residue	50.0	40.0 to 60.0		

4.	TEST MEASUREMENT:						
	a) test initiated immediately after preparing and trimming sample						
b) sample conditioned by applying							
i) a recommended 10 cycles at 10 rad/s							
ii) a required strain 8 to 16 cycles at 10 rad/s							
c) test measurement acquired by repeating the:							
i) recommemded strain for 10 more cycles							
ii) required strain for 8 to 16 more cycles							
	c) data obtained for a second of cycles reduced to produce a value for complex						
	modulus and phase angle						
	d) when testing at more than one frequency testing started at the lowest temperature						
	e) all testing completed within 4 hours						
	<u>REPORT</u>						
	SAMPLE IDENTIFICATION: sample name, type, code number, source etc						
2.	INSTRUMENT USED:						
	a) whether at constant stress or strain rheometer						
	b) type of environmental chamber						
	c) report observed modulus and whether or not strain levels were outside the linear						
	range						
	d) report modulus at recommended stress or strain levels						
	e) report if data was obtained oudside the linear region						
	TEST PLATE DIAMETER: to nearest 0.1 mm						
	TEST GAP: to nearest 1 µm						
	TEST TEMPERATURE: to nearest 0.1°C						
	TEST FREQUENCY: to nearest 0.1 rad/s						
	STRAIN AMPLITUDE OR TORQUE: to nearest 0.01 % or nearest mN.m respectively						
9.	COMPLEX MODULUS: (G*) for 10 measurements, kPa to three significant figures						
10	PHASE ANGLE: phase angle (δ) for second 10 cycles, report to nearest 0.1°C						
11	QUIPMENT MANUFACTURER:						
12	MODEL:						

REMARKS:			